


Microscope Safety

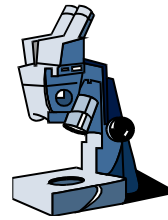

1. Always carry the microscope with 1 hand under the _____ and 1 hand on the _____.
2. _____ the cover, _____ the cord, and plug it into the _____ outlet.
3. The _____ position for the microscope is always stage _____ and the _____ power objective over the stage.
4. Place the _____ slide on the stage, fasten the _____, raise the stage with the _____ adjustment knob, and adjust the _____ by rotating the wheel to let in the best amount of _____.
5. 1st View - adjust with the _____ adjustment knob and rotate the eyepiece _____ so that it points to the clearest specimen.
6. 2nd View - lower the _____, rotate the revolving _____ to the _____ power objective, raise the _____, and adjust with the _____ adjustment knob (also, adjust the _____ pointer as needed).
7. 3rd View - lower the _____, rotate the _____ nosepiece to the _____ power objective, _____ and _____ raise the stage, and adjust with the _____ adjustment knob (also, adjust the eyepiece pointer as needed).
8. Complete all of the laboratory _____ including _____ work.
9. Return the microscope to the _____ position and place the _____ slide in the appropriate _____.